OneLab How to Set Up a Meta Quest Headset Follow these steps to set up your headset: Start with the comprehensive guide from Meta 🔿 Meta Log in or create a Meta account 1. Create a new Meta account at Continue with Facebook auth.meta.com O Continue with Instagram and select the "Continue Continue with email with Email" button. Android 2. Install Meta app on GET IT ON your mobile device. Google Play Search for Meta App in your mobile device app store. Once installed, sign in with the **iPhone** Note: account you created using your email. d on th Follow the prompts. You will need a Wi-Fi App Store connection and a mobile device (phone or tablet) to complete this process. 6. Adjust 3. Charge your head straps headset and distance with the included USB-C between lenses cord; be sure each of the controllers has AA batteries installed 4. Power on Headset 5. Choose Press and hold power your preferred button until you hear a chime language (Meta Quest default language is English)

Questions? Email VR@cdc.gov or visit reach.cdc.gov/onelabvr

How to Set Up a Meta Quest Headset



8. Ack

8. Acknowledge license agreement

OneLab

9. Headset software will update

10. Connect headset to mobile Meta App

Open the Meta app and select the Menu button at the bottom, then select the Devices button and follow the prompts for your Meta Quest headset. If your headset does not connect automatically, you may need to enter the PIN displayed in the headset.

11. Set up 'Guardian' boundary

Follow the prompts to set the floor position and define a boundary. Give yourself at least 6.5 square feet of open space. You also have a stationary option if you have limited space.



New to VR?

Try Meta's First Steps app to get some practice using the controllers, navigation, and other tips. To find new apps go to the Meta Store.

Bluetooth

For Tech Support

please check Meta's support at **meta.com/help/**